



www.mypols.de  
support@mypols.de  
T +49(0)7531 122 965 00

## Taq DNA polymerase

#1101 (hotstart), Manual Version 2, January 25, 2021

Store at -20°C

### Contents

Taq DNA polymerase is supplied as a 5 U/μl solution containing glycerol. It comes together with a 10x optimized reaction buffer.

### Description

Taq DNA polymerase is supplied together with the 10x Taq reaction buffer.

The reaction buffer has been specifically designed for optimal PCR performance and polymerase activity. Taq DNA polymerase can also be used for realtime cycling, when adding a suitable realtime dye or a fluorescent probe.

### Applications

- Standard PCR
- Realtime-PCR (addition of suitable dye required)
- Primer extension reactions
- TA cloning
- 3'A-tailing of blunt ends
- Screening / High-throughput PCRs

### Troubleshooting

How can I optimize the PCR reaction conditions?

1. The annealing temperature can usually be optimized. Try a temperature gradient and determine the best annealing temperature, which yields in the cleanest product.
2. Add a gradual amount of betaine 0-1M or DMSO 0-7.5% to the reaction mix and select for the cleanest product and the highest yield.
3. Try to shorten the extension and annealing time. Too long and too many cycles may lead to over-amplification and side-products.

### References

Isolation, characterization, and expression in *Escherichia coli* of the DNA polymerase gene from *Thermus aquaticus*. *J. Biol. Chem.* 1989; 264 (11):6427-6437. F. C. Lawyer, S. Stoffel, R. K. Saiki, K. Myambo, R. Drummond, and D. H. Gelfand.

Fidelity of DNA synthesis by the *Thermus aquaticus* DNA polymerase. *Biochemistry* 1988; 27(16): 6008-6013. K. R. Tindall, T. A. Kunkel.

Primer-directed enzymatic amplification of DNA with a thermostable DNA polymerase. *Science* 1988; 239(4839): 487-491. R. K. Saiki, D. H. Gelfand, S. Stoffel, S. J. Scharf, R. Higuchi, G. T. Horn, K. B. Mullis and H. A. Erlich.

## Recommendations for PCR/ Reaction Setup

### PCR Mix

Component	Volume	Final concentration
Primer forward (10 μM)*	1 μl	0.2 μM (0.05-1 μM)
Primer reverse (10 μM)*	1 μl	0.2 μM (0.05-1 μM)
dNTPs (2 mM)	5 μl	200 μM
Buffer (10x)	5 μl	1x
Taq DNA polymerase 5 U/μl	0.25 μl	1.25 U/reaction
Template/Sample extract	x μl	<1000 ng** DNA
Nuclease-free water		up to 50 μl total vol.

\* Primers should ideally have a GC content of 40-60% typically

\*\*Suggested template concentration should be about 1 ng - 1 μg (genomic DNA) or 1 ng - 1 pg (plasmid/viral DNA).

### Typical 3-step PCR protocol

Initial denaturation	95°C	2 min	} 25-40 cycles
Denaturation	95°C	15 sec	
Annealing*	54-72°C	30 sec	
Extension	72°C	1 min/1000 bp	
Hold	<10°C		

\* Typically, the annealing temperature is about 3-5°C below the calculated melting temperature of the primers used.

### Important notes

- Keep all components on ice.
- Spin down and mix all solutions carefully before use.
- Primers should ideally have a GC content of 40-60%.
- Suggested template concentration should be about 1 ng - 1 pg (plasmid/viral DNA) or 1 ng - 1 μg (genomic DNA).



### Quality Control Assays

Taq DNA polymerase is tested for successful PCR performance. A 92 bp fragment (beta-actin gene) was amplified from human genomic DNA and analysed by agarose gel electrophoresis. The activity of Taq DNA polymerase has been monitored and adjusted to a specific DNA polymerase activity using an artificial DNA template and DNA primer. Enzyme-concentration has been determined by protein-specific staining. Please inquire more information at info@mypols.de for the lot-specific concentration. No contamination has been detected in standard test reactions.

### Material Safety Data (MSDS)

According to OSHA 29CFR1910.1200, Australia [NOHSC:1005, 1008 (1999)] and the EU Directives 67/548/EC, 1999/45/EC and 1272/2008 (CLP Regulation) any products which do not contain more than 1% of a component classified as dangerous or hazardous nor more than 0.1% of a component classified as carcinogenic, do not require a MSDS. However, we recommend the use of gloves, lab coats and eye protection when working with these or any other chemical reagents. myPOLS Biotec takes no liability for damage resulting from handling or contact with this product. This product is not hazardous, not toxic, not IATA-restricted. Product is not from human, animal or plant origin. The source of the product is recombinant protein expression in *E. coli*. The product is for research use only and may be used for *in-vitro* experiments only.